



Gravity Well Instruction Manual

© Magenta LLC. 2021

Using the Vessels

Vessels can be used like a Magenta box or other reusable vessel with semi-solid or with liquid medium.

Always autoclave with the lids on and seal the edges with 3M Micropore tape or similar sealing film.

Using the Gravity well

Step 1

Place a strip of regular printer paper along the bottom of the vessel. Cut the paper to 65mm x 210mm or (2.5" x 8.25").

Step 2

Place the Gravity Well on one side of the vessel.

Step 3

Fill with 100mL of liquid media.




Step 4

Autoclave, then let cool.

Step 5

Tip the Gravity Well to refill the reservoir.

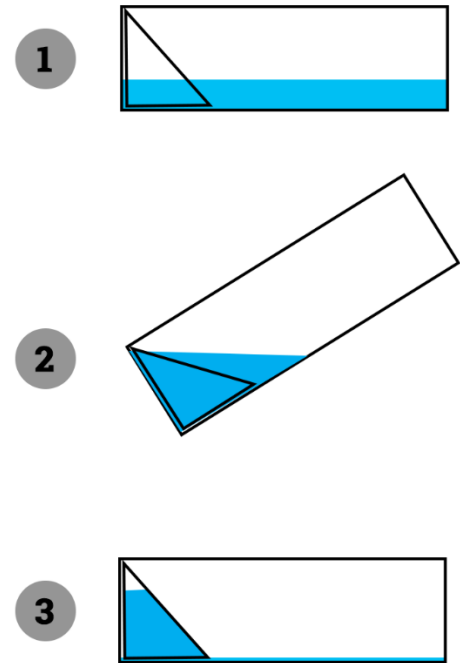

Step 6

Place shoots along the paper on the bottom of the vessel.



Caution: Vibration and motion can cause the liquid layer to become too thick. If this occurs, you can tip the vessel to refill the reservoir again, but go slowly so plants don't enter the reservoir.

Always use the Gravity Well with Ventilated lids.

Troubleshooting

If you are experiencing any problems, please use the contact form at www.wevitro.com