



Photoblaster

Instruction Manual

© Magenta LLC. 2021

Using the Vessels


Vessels can be used like a Magenta box or other reusable vessel with semi-solid or with liquid medium.

Always autoclave with the lids on and seal the edges with 3M Micropore tape or similar sealing film.

Assembling the LED Lights

To assemble the Photoblasters, first remove the LED light bar from the plastic packaging.

Remove the beige plastic LED clips from the vessels.

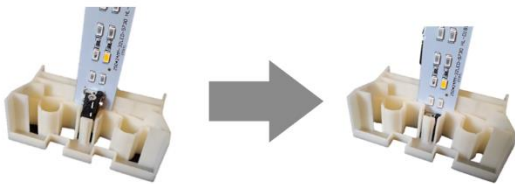

Next, place the LED clips on a flat surface. Push the LED light bar into the center slot of the LED clips. You will hear a click when the LED bars have snapped into place. Some force may be required. Repeat this step for the other end.

Connecting the Photoblasters



Plug the included AC power adapter into the wall.



Then plug in the LED dimmer module onto the end of the AC power adapter.



Connect the LED dimmer to the first Photoblaster. Use the included double-ended cables to connect Photoblasters together.

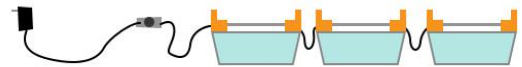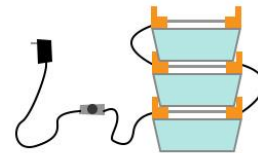

Only use 5 Photoblasters in a single chain.

Troubleshooting

If you are experiencing any problems, please use the contact form at www.wevitro.com