


Best Practices and Care Sheet

- Always autoclave vessels with the lid on to avoid warping
- Autoclave at 121°C at 118 kPa for 20-40 minutes
- Seal vessels with 3M Micropore tape (after adding plants)
- To remove lid easily, grasp one end and wiggle:

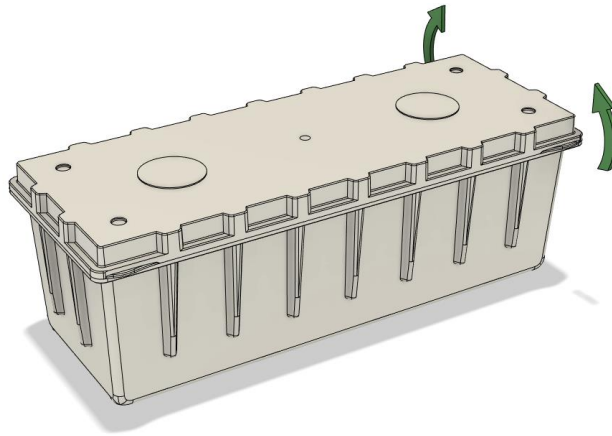

To avoid clouding / scabbing of the plastic

- Avoid glassware detergents (Sparkleen and alkaline detergents)
– use plastic-safe detergents for cleaning*
- Do not use heated dry cycle on dishwashers – hand dry with a dish towel or -
- Use polycarbonate-safe rinse agent or a vinegar rinse

*: L900 Nalgene, Joy, Palmolive liquid, Top Job